

TwistAmp® Liquid exo/exo RT

Quick Guide

Part number: LQEXOQG01

Revision 03V4

TwistAmp® Liquid exo/exo RT Quick Guide

Part number: LQEXOQG01 / Revision 03V4

Please see **instruction** and **assay design manuals** at twistdx.co.uk for information regarding components and storage, assay design, and detailed use.

Instructions are based on 50 µl reaction volumes; if using a different volume, quantities should be adjusted appropriately.

Primer screen set-up (single-plex)¹

1. Add 2.1 µl of each primer and 0.6 µl of exo probe at 10µM concentration to 0.2 ml PCR tubes.

2. Prepare a pre-master mix (per reaction) in the order below:

2x Reaction Buffer	25 µl
dNTPs ² + water to	8.2 µl
(7.2 µl if adding RT ^{3,4})	
10x Probe E-mix	5 µl

Vortex and spin briefly.

3. To the pre-master mix, add 2.5 µl 20x Core Reaction Mix⁵ (per reaction) to tube lid. Mix by 10x full inversions and spin briefly. If using RNA, add 1 µl 50x RT (per reaction) to tube lid.

4. Add 1 µl 50x Exo (per reaction) to tube lid. Mix by 10x full inversions and spin briefly. Master mix is now complete⁶. Pipette mix before use.

5. Add 41.7 µl^{3,4} master mix to primers and probe prepared in tubes (step 1) and pipette mix.

6. Add 2.5 µl of 280mM MgOAc (supplied) and 1 µl template to tube lids^{3,4}. DNA/RNA and MgOAc should be kept separate in the tube lid prior to spin-down. Spin in MgOAc/ template and mix well (6x inversions) to start reaction. Spin briefly.

Warning: RPA reactions start as soon as MgOAc is added.

7. Place reactions in a fluorometer and start run: 37-42°C, 20 minutes. For low template copies, remove strip after 4 mins (5 mins for RNA), mix by 6x full inversions and spin briefly, replace in fluorometer.

TwistAmp® Liquid exo/exo RT Quick Guide

Part number: LQEXOQG01 / Revision 03V4

Template screen set-up (single-plex)¹

1. Prepare a primer and probe pre-master mix (per reaction) in the order below:

2x Reaction Buffer	25 µl
dNTPs ² + water to	8.2 µl
(7.2 µl if using RT ^{3,4})	
10x Probe E-mix	5 µl
Primer A (10µM)	2.1 µl
Primer B (10µM)	2.1 µl
Probe (10µM)	0.6 µl

Vortex and spin briefly.

2. Add 2.5 µl 20x Core Reaction Mix⁵ (per reaction) to tube lid. Mix by 10x full inversions and spin briefly.

3. If using RNA, add 1 µl 50x RT (per reaction) to tube lid. Add 1 µl 50x Exo (per reaction) to tube lid. Mix by 10x full inversions and spin briefly. Master mix is now complete⁶. Pipette mix before use.

4. Add 46.5 µl^{3,4} master mix to 0.2 ml PCR tubes.

5. Add 2.5 µl of 280mM MgOAc and 1µl template to tube lid^{3,4}. DNA/RNA and MgOAc should be kept separate in tube lids prior to spin-down.

Spin in MgOAc/template and mix well (6x inversions) to start reaction. Spin briefly.

Warning: RPA reactions start as soon as MgOAc is added.

6. Place reactions in fluorometer and start run: 37-42°C, 20 minutes. For low template copies, remove strip after 4 mins (5 mins for RNA), mix by 6x full inversions and spin briefly, replace in fluorometer.

- ¹ See manual for multiplexing.
- ² Suggested final concentration of 1.8mM (total) dNTPs. Optimisation is recommended.
- ³ If amplifying RNA, users may wish to add RNase Inhibitor together with the template and adjust volumes accordingly.
- ⁴ Volumes should be adjusted if adding more/less template and/or MgOAc.
- ⁵ Warm to room temperature and pipette mix slowly to ensure homogeneity.
- ⁶ Master mix may appear cloudy, this is normal.

Notice to purchaser

Licence, Use Restrictions and Limitations of Liability

Definitions. As used in this section, “kit” means the items described in this manual (the “Manual”) and supplied by TwistDx to a purchaser (the “Recipient”). “Materials” means all biological and chemical materials supplied as part of the kit. “Information” means all written information supplied as part of the kit, information relating to the kit made available through TwistDx’s website, and any verbal or written information concerning the kit or its use provided by any employee or agent of TwistDx.

Limitations on Use and Distribution of the Material and Information

Recipient acknowledges and agrees that the Materials and Information are proprietary to TwistDx, may be covered by claims of patents or patent applications owned by TwistDx or its affiliates and are supplied subject to the following restrictions: The Materials and Information are non-exclusively licensed to Recipient solely for non-commercial internal research purposes and for applications other than the sequential determination of the identity and relative order of at least 200,000 total nucleotides in a single run on a sequencing apparatus. Any *in vitro* diagnostic use of the Materials and Information or any use for diagnosing or monitoring any medical condition in a human is expressly prohibited.

No Warranty; Limitation of Liability

Recipient understands and agrees that the materials are experimental in nature and that the Materials and Information are provided without any warranty as to results, merchantability, fitness for a particular purpose or non-infringement of any patent or other intellectual property right, and without any other representation, warranty or condition, express or implied. TwistDx shall not be liable in connection with the Materials, Information or any breach of this agreement under any contract, negligence, strict liability or other theory for (a) loss of revenues, loss of profits, or loss or inaccuracy of data, including test results, regardless of how such damages are characterised (b) for the cost (including procurement costs) of substitute goods, services or technology, or (c) for any special, indirect, incidental or consequential damages. In no event will TwistDx’s aggregate liability under this agreement exceed one hundred dollars (US\$100). Recipient understands that its use of the materials and/or the information in connection with its activities are entirely at its risk.