## Notice to purchaser

Absent a written agreement between you and TwistDx Limited to the contrary, the supply of this manual ("Manual") and the product described herein ("Product") is subject to the TwistDx Limited Terms and Conditions of Supply available at https://www.twistdx.co.uk/terms-of-supply, including, without limitation, the use restrictions, warranty disclaimer and limitation of liability set forth therein.

#### **Use Restrictions**

FOR RESEARCH USE ONLY. NOT FOR USE IN DIAGNOSTIC PROCEDURES.

The Products are for non-commercial internal research use only. The Products must not be used for any *in vivo* testing, nor with humans or animals for any clinical diagnostic purposes. See the TwistDx Limited Terms and Conditions of Supply available at https://www.twistdx.co.uk/terms-of-supply for additional use restrictions.

#### No Warranty

TwistDx Limited makes no warranty with respect to the Product and excludes any warranties implied by law to the fullest extent permitted by law. THE PRODUCT IS PROVIDED WITHOUT ANY WARRANTY AS TO RESULTS, MERCHANTABILITY, FITNESS FOR A PARTICULAR PURPOSE OR NON-INFRINGEMENT OF ANY PATENT OR OTHER INTELLECTUAL PROPERTY RIGHT, AND WITHOUT ANY OTHER REPRESENTATION, WARRANTY OR CONDITION, EXPRESS OR IMPLIED, EXCEPT AS STATED IN THE PRECEDING SENTENCE. See the TwistDx Limited Terms and Conditions of Supply available at https://www.twistdx.co.uk/terms-of-supply for additional terms and conditions.



## TwistAmp<sup>™</sup> nfo Kit

# Quick Guide

Part number: INTANFO Revision 3



Abbott House, Vanwall Business Park, Vanwall Road, Maidenhead, SL6 4XE, UK E info@twistdx.co.uk | twistdx.co.uk

> © 2019 Abbott. All rights reserved. All trademarks referenced are trademarks of either the Abbott group of companies or their respective owners.

Please see instruction and assay design manuals at twistdx.co.uk for information regarding kit components and storage, assay design, detailed use and multiplexing.

### Set-up (single-plex)

1. Prepare reaction mix in 1.5ml tube:

Primer A (10µM)	2.1 µl
Primer B (10µM)	2.1 µl
TwistAmp <sup>®</sup> nfo probe (10µM)	0.6 µl
Primer Free Rehydration buffer	29.5 µ
Template and water to	13.2 µ
(Total volume	47.5 μ
Vortex and spin briefly.	

- 2. Add reaction mix to a TwistAmp® nfo reaction. Pipette to mix.
- Add 2.5 µl of 280mM Magnesium Acetate (MgOAc) (supplied) and mix well to start reaction.

### **Note:** RPA reactions start as soon as MgOAc is added.

- Incubate at 40°C for 20 minutes. For low template copy number, remove strip after 4 minutes, vortex and spin briefly, replace in heating device.
- After 20 minutes, for analysis by lateral flow, dilute reaction products as specified by the lateral flow consumable of choice (see instruction manual).

**Note:** If tubes are opened after amplification there is a great risk of contamination of work surfaces with amplicon. Ensure that appropriate avoidance measures are taken.

TwistAmp® exo probes can also be used with this kit (see instruction manual).

### Kit Positive Control Set-up (single-plex)

- 1.Prepare reaction mix in 1.5ml tube:Positive control primer mix8 μlPrimer Free Rehydration buffer29.5 μlPositive control DNA template1 μlWater9 μl(Total volume47.5 μl)Vortex and spin briefly.
- 2. Add reaction mix to a TwistAmp® nfo reaction. Pipette to mix.
- Add 2.5 µl of 280mM Magnesium Acetate (MgOAc) (supplied) and mix well to start reaction.

### Note: RPA reactions start as soon as MgOAc is added.

- Incubate at 39°C for 20 minutes. For low template copy number, remove strip after 4 minutes, vortex and spin briefly, replace in heating device.
- After 20 minutes, for analysis by lateral flow, dilute reaction product 1/50 with PBST and load 10 µl onto the sample pad of Milenia HybriDetect 1 strip.
- 6. Place strip in PBST running buffer.

**Note:** If tubes are opened after amplification there is a great risk of contamination of work surfaces with amplicon. Ensure that appropriate avoidance measures are taken.

TwistDX	PMS 185 C	Part Number: INTANFO
nfo Kit QG	Red	Revision 3
Size:	PMS 7541 C	Date of Revision:
8.268 in x 5.827 in	Gray 1 - 10%	3.1 2019/01